



# Laser Controller

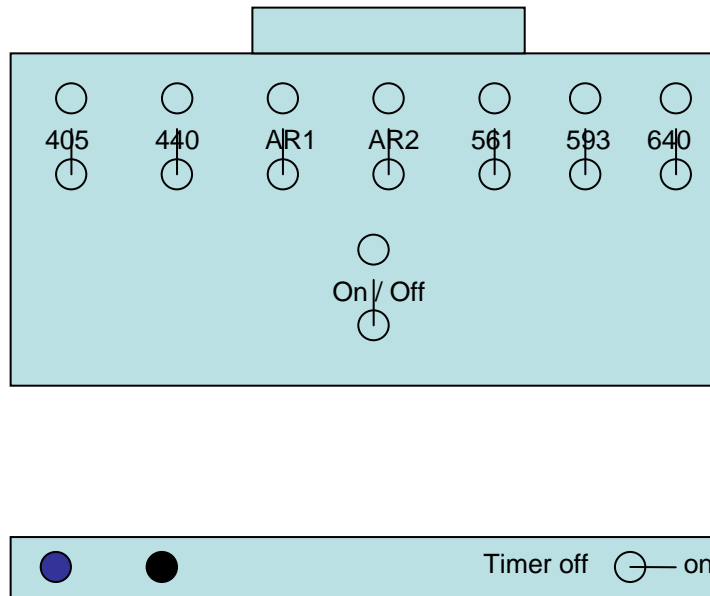

1. Select the laser you want to use
2. Toggle the on/off switch to on
3. For use without timer check that the timer is switched to the off position

# Timer Box

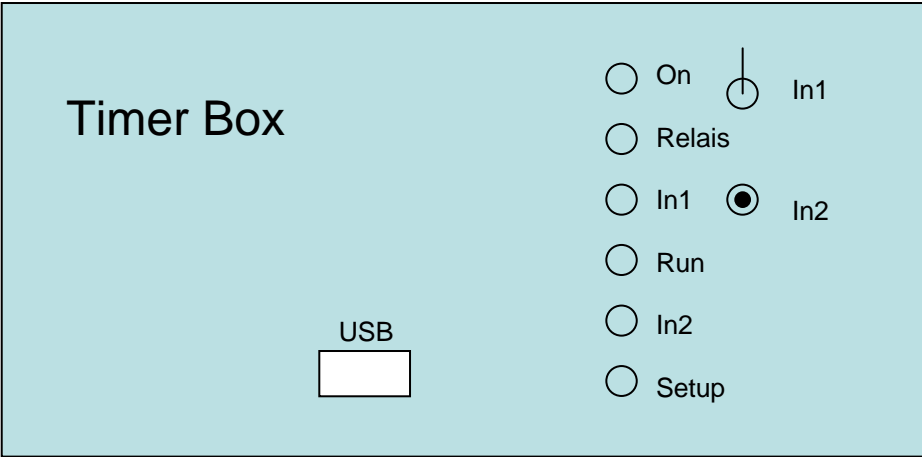

## Setup timer

1. Plug-in USB cable.
2. Start HB626 software (fig. 1).
3. Select Fn5 (fig. 2) en fill in time in seconds.
4. Select “Einstellung abschicken” and Figure 3 should appear.
5. Select OK and close program.
6. Remove USB cable. The timer starts immediately. To reset the starting time you can switch In1 off and on again.

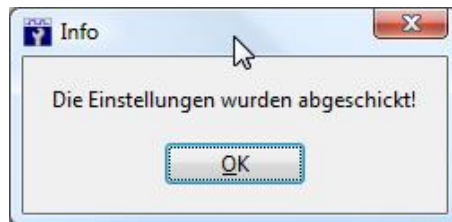


Figure 3

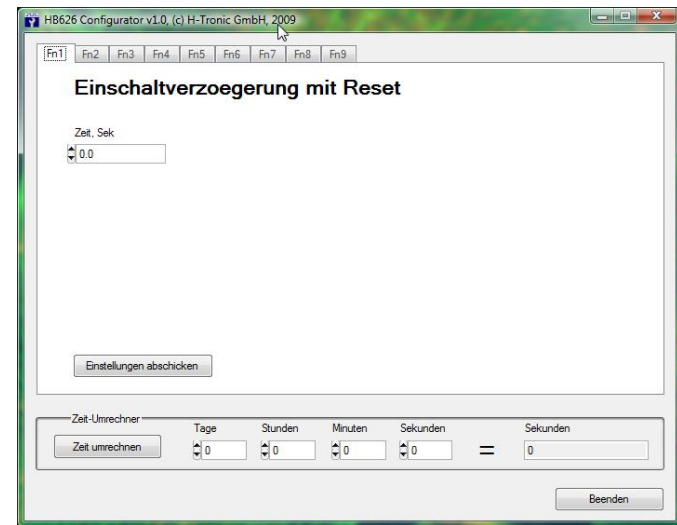


Figure 1

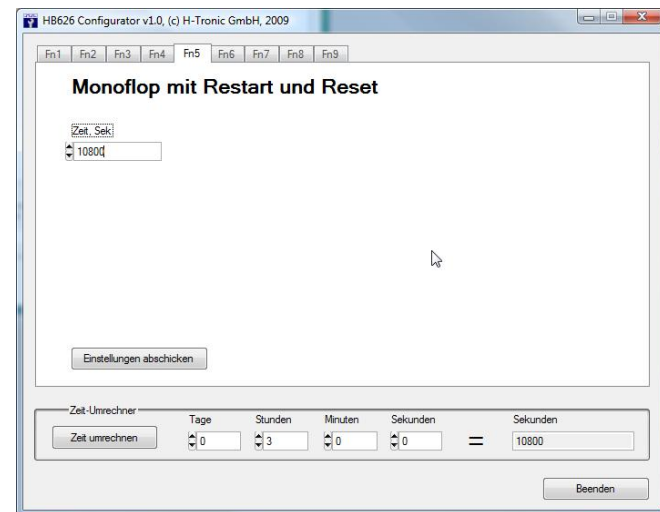


Figure 2

7. To use the timer for the lasers toggle the timerswitch on the laser controller to on.